

EZ PANGASIUS

A Fast, Sensitive and Simple Test for the Detection of Cooked or Raw Pangasius Filets

- + QUICKLY DETECTS PANGASIUS SPECIES
- + SIMPLE - SMALL- PORTABLE
- + TEST RAW AND COOKED PRODUCTS
- + TEST WHOLE FILETS
- + FOR FIELD USE
- + EASY TO USE

With the EZ Pangasius™ Test, Measure out one level spoonful of the diced filet using the spoon provided, or weigh out 0.5g of the sample on a scale

1.



then add the sample to the pangasius extraction solution and mix.

2.



Place a few drops of the sample extract into a test tube.

3.




The EZ Pangasius™ test strip is placed into the test tube to absorb the sample extract.

4.



After 10 minutes, the test strip can be read for the presence of pangasius.

5.

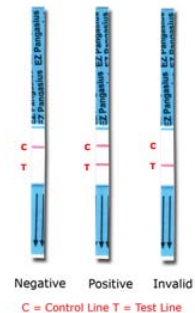

ELISA Technologies, Inc.



EZ Pangasius™ - 10 Test Kit 510EZP

Call us Toll-Free
1-866-222-4729
For Pricing and Shipping